# VWR MiniArray

## INSTRUCTION MANUAL

### European Catalogue Number(s):

**Euro Plug: VWRI521-2849**

**UK Plug: VWRI521-2850**

**Swiss Plug: VWRI521-2851**

Version: 3

Issued: 2011-09-15

## Legal Address of Manufacturer

### **Europe**

VWR International bvba  
Researchpark Haasrode 2020  
Geldenaaksebaan 464  
B-3001 Leuven  
+ 32 16 385011  
<http://be.vwr.com>

Country of origin    South Korea

# Table of Contents

<b>Safety information</b>	<b>2</b>
<b>Package Contents</b>	<b>2</b>
<b>Intended Use</b>	<b>2</b>
<b>Product specifications</b>	<b>2</b>
<b>General Operating Instructions</b>	<b>3</b>
<b>Maintenance</b>	<b>3</b>
<b>Questions and service</b>	<b>3</b>
<b>Technical service</b>	<b>4</b>
<b>Warranty</b>	<b>4</b>
<b>Equipment Disposal</b>	<b>5</b>

## Safety Information

**NEVER** use the centrifuge in any manner not specified in these instructions

**NEVER** operate the centrifuge without a rotor properly attached to the shaft.

**NEVER** put hands in the rotor area unless the rotor is completely stopped

**NEVER** move the centrifuge while the rotor is spinning

**NEVER** use solvents or flammables near this or other electrical equipment

**NEVER** centrifuge flammable, explosive, or corrosive materials

**NEVER** centrifuge hazardous materials outside of a hood or proper containment facility

**NEVER** spin slides that do not fit easily into the cassettes.

**NEVER** autoclave. Autoclaving can cause the polycarbonate to lose its strength over time.

**ALWAYS** load the rotor symmetrically. Each tube should be counter balanced by another tube. Weight of individual tubes may not exceed 3.0 g.

**ALWAYS** locate the centrifuge within easy access to an electrical outlet

**ALWAYS** load the rotor symmetrically. Each slide should be counterbalanced by the second slide.

## Package Contents

Description	Quantity
Mini Centrifuge with 2 place slide rotor	1
Power cord	1
Slide holders	2

## Intended use

*The MiniArray is a personal slide drying centrifuge. This centrifuge will dry 2 slides without splashing. The unit is supplied complete with the 2 place rotor and 2 reusable cassettes.*

*Additional cassettes may be purchased separately. Please read these operating instructions thoroughly before attempting to operate the centrifuge.*

## Product Specifications

<b>Dimensions</b>	132 X 155 X 110 mm
<b>Electrical</b>	230V, 50/60Hz
<b>Fuse</b>	0.5A, 250 V (F)
<b>Max Speed</b>	4,000 rpm
<b>Max RCF</b>	500-850 x g
<b>Power Consumption</b>	15W

## General Operating Instructions

Insert a slide into each of the cassettes on the rotor. **Note: Always run 2 slides at once.** Use a blank slide as a balance if only one slide needs to be dried. To begin a run, simply close the lid of the centrifuge. No other controls are required. Typically a 10 second spin is required to completely dry the microarray slides. To stop rotation, press down on the lid release tab on the front of the unit or put the switch on the side to the "off" position. Pressing the tab will release the lid lock and the rotor will slow gradually to a complete stop. When the rotor has come to a complete stop, open the lid and remove the slides. The cassettes can be removed from the rotor for easy disposal of the liquid that collects at the bottom of the cassettes.

**Caution:** Do not attempt to open the lid or remove samples until the unit has come to a complete stop. After the rotor has stopped, the lid may be opened by grabbing it with the thumb on the front and fingers on the back then lifting the lid back on the hinge.

Safe operation of the centrifuge requires that the rotor be loaded in a balanced fashion. Failure to load the rotor properly presents a hazardous condition and may damage the centrifuge. Always run 2 slides for proper rotor balance.

## Maintenance

Before using any cleaning or decontamination methods except those recommended by the manufacturer, check with the manufacturer that the proposed method will not damage the equipment.

To clean the centrifuge: use a damp cloth and mild, noncorrosive detergent (pH<8). After cleaning, ensure all parts are dried thoroughly before attempting to operate the unit. Do not immerse the centrifuge in liquid or pour liquids over it.

The recommended cleaning method for the cassettes is with a mild non-alkaline detergent. For decontamination of the internal part of the cassettes the manufacturer recommends using Ultrasonic Cleaning Units (it is very effective for destroying DNA – autoclaving doesn't effectively destroy DNA).

**The cassettes are not resistant to alkaline detergents, or acids**

## Questions and Service

Should you have a question about the centrifuge or require service for the unit, please contact your VWR International representative. Do not send in a unit for service without first calling to obtain a repair authorization number and a decontamination form. Should the centrifuge require service, be sure to decontaminate it and enclose the decontamination form. The unit should be properly packed to avoid damage. Any damage resulting from improper packaging shall be the responsibility of the user.

## Technical service

### Web Resources

Visit the VWR's website at [www.vwr.com](http://www.vwr.com) for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

**Contact us** For information or technical assistance contact your local VWR representative or visit.

[www.vwr.com](http://www.vwr.com).

## Warranty

**VWR International** warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of purchase. If a defect is present, VWR will, at its option, repair, replace, or refund the purchase price of this product at no charge to you, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear.

For your protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

## Equipment disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

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Typically a 10 second spin is required to completely dry the microarray slides. To stop rotation, press down on the lid release tab on the front of the unit or put the switch on the side to the "off" position. Pressing the tab will release the lid lock and the rotor will slow gradually to a complete stop. When the rotor has come to a complete stop, open the lid and remove the slides. The cassettes can be removed from the rotor for easy disposal of the liquid that collects at the bottom of the cassettes.

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